



**AZ Instrument Corp.**

Web site: <http://www.az-instrument.com.tw>  
E-mail: [info@az-instrument.com.tw](mailto:info@az-instrument.com.tw)

Tel: 886-4-25326668 (rep.)  
Fax: 886-4-25326593

---

Edit by Naima 11 Feb. 2020

## **Question**

**What to do when display shows pH 7 at every calibration points?**

先問序號，若查序號發現已出貨超過一年則回

## **Reply:**

After look up on the serial number, due to the item already been shipped out for more than one year, we assume that the reason is because of electrode aging.

Nevertheless, you can still try to clean the probe by immersing the electrode in tap water for 10-15 minutes, then rinse thoroughly with distilled water, to see if it's still available.