



Edit by Naima 30th Sep. 2020

What's the procedure of pH calibration? When should I re-calibrate my pH meter?

AZ86851 pH pen offer 3 points calibration available, please always do pH 7 calibration first, and then pH 4 or pH 10 depends on the measuring solution is closer to which range for better accuracy

And in the following cases, the meter must be re-calibrated before use:

1. The electrode hasn't been used for a long time or the electrode is brand new.
2. After measuring strong acid (pH<2) or strong base (pH>12) solutions.
3. After measuring fluoride-containing solution and strong organic solution
4. There is a big difference between the temperature of the test sample and the temperature of the buffer solution that is used in the last calibration.